

Contents

1. Description
 - 1.1 Background information
 - 1.2 Applications
2. Protocol
 - 2.1 Preparation
 - 2.2 Instructions for use

1. Description

Components	100 mL MACS® Tissue Storage Solution
Sterility	Bioburden tested.
Storage	Store protected from light at 2–8 °C. Do not freeze. The expiration date is indicated on the label.

1.1 Background information

The storage of organs or tissue samples may be required if material has to be collected from multiple origins, at different time points, or if it can not be directly processed. During this storage period necrosis and apoptosis must be avoided and cells like stem cells need to be prevented to lose their functional capacities. Immune cells may be activated causing a bias in downstream analysis during storage. The MACS® Tissue Storage Solution has been developed to allow an optimized storage of fresh organ and tissue samples without showing background effects like cell activation or apoptosis induction. It has been tested and validated on a variety of human and murine tissues including tumors, skin, heart, spleen, brain, and skeletal muscle.

1.2 Applications

- Storage of fresh organ and tissue samples
- Shipping of fresh organ and tissue samples

2. Protocol

2.1 Preparation

- ▲ The MACS Tissue Storage Solution is formulated without antibiotic or antifungal components. If required, supplement with these ingredients at standard cell culture concentrations prior to use.
- ▲ All procedures should be carried out using sterile techniques in a biosafety cabinet.
- ▲ Pre-chill MACS Tissue Storage Solution to <8 °C before use.

2.2 Instructions for use

1. Transfer sample directly into the pre-chilled MACS Tissue Storage Solution (<8 °C). If the sample has already been placed in other buffers or media, replace the complete liquid with the pre-chilled MACS Tissue Storage Solution. The sample must be completely covered with the MACS Tissue Storage Solution.
2. (Optional) For shipping purpose, transfer sample into the MACS Tissue Storage Solution in a closable tube. Avoid air bubbles to prevent air contact for prolonged time periods.
3. Store or ship sample at 2–8 °C.
4. At the end of storage, remove sample from the MACS Tissue Storage Solution and process directly with downstream application, e.g., dissociation into single-cell suspensions by using the gentleMACS™ Dissociator or gentleMACS Octo Dissociator.

All protocols and data sheets are available at www.miltenyibiotec.com.

Warranty

The products sold hereunder are warranted only to be free from defects in workmanship and material at the time of delivery to the customer. Miltenyi Biotec GmbH makes no warranty or representation, either expressed or implied, with respect to the fitness of a product for a particular purpose. There are no warranties, expressed or implied, which extend beyond the technical specifications of the products. Miltenyi Biotec GmbH's liability is limited to either replacement of the products or refund of the purchase price. Miltenyi Biotec GmbH is not liable for any property damage, personal injury or economic loss caused by the product.

MACS is a registered trademark and gentleMACS is a trademark of Miltenyi Biotec GmbH.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research use only and not for diagnostic or therapeutic use.

Copyright © 2013 Miltenyi Biotec GmbH. All rights reserved.